
Bench Top Extraction Arms

FX₂ Original/Chem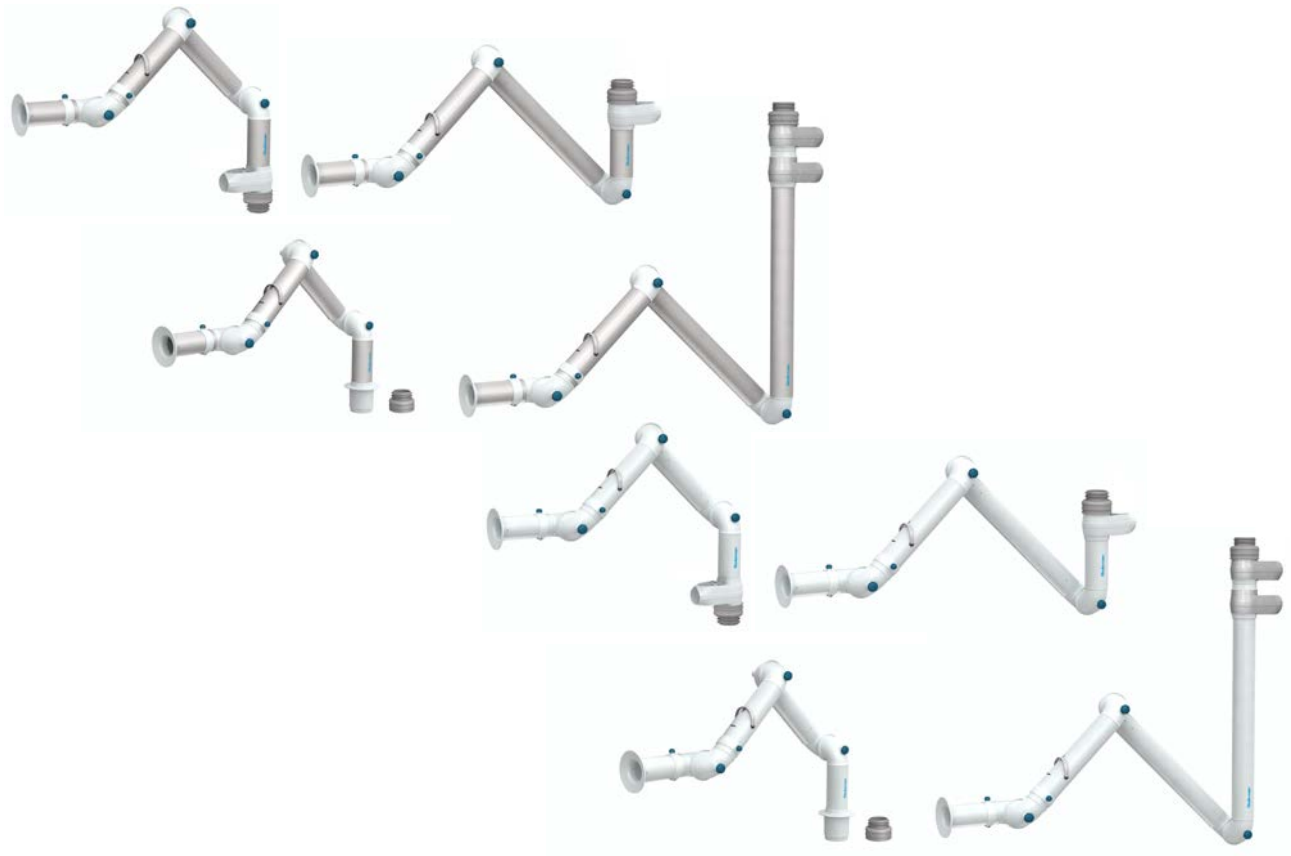

Original instruction manual

EN INSTRUCTION MANUAL

Translation of original instruction manual

CS NÁVOD K OBSLUZE
DE BEDIENUNGSANLEITUNG
ES MANUAL DE INSTRUCCIONES
FI KÄYTTÖOHJEET
FR MANUEL D'INSTRUCTION

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EN English

Declaration of conformity

We, AB Ph. Nederman & Co., declare under our sole responsibility that the Nederman product:

FX₂ Original/Chem (Part No. **, and stated versions of **) to which this declaration relates, is in conformity with all the relevant provisions of the following directives and standards:

Directives

2006/42/EC.

Standards

EN ISO 12100.

The name and signature at the end of this document, is the person responsible for both the declaration of conformity and the technical file.

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Chem: 70603144-70603844, 70605644-70606444

AB Ph. Nederman & Co.
P.O. Box 602
SE-251 06 Helsingborg
Sweden




Dorinel Lapadat
Manager Product Centre
2017-03-17

Figures

1

Duct Connectors

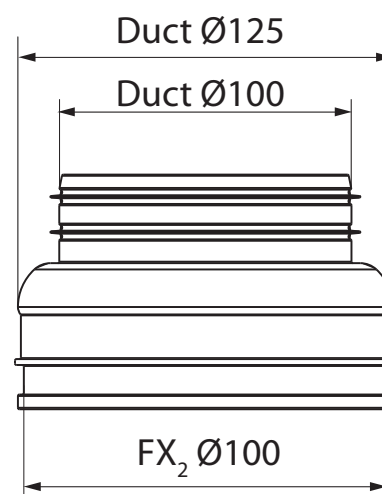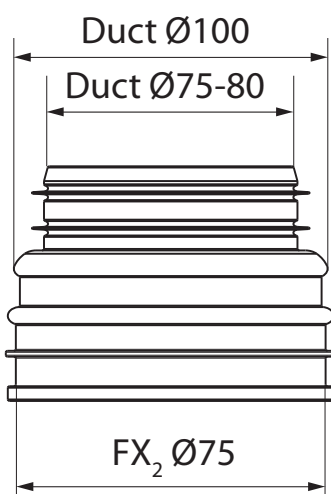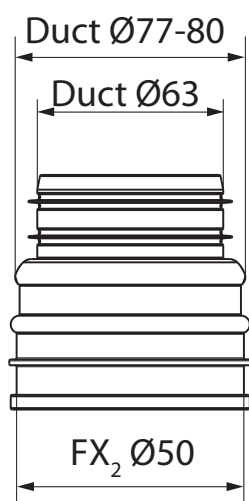2

Technical Data

Ø50 mm		
FX ₂ – D50 – Up/Down - Low	FX ₂ – D50 – Up - PLUS	FX ₂ – D50 – Up/Down - Through
1,9 – 2,8 kg	2,5 – 3,4 kg	1,9 – 2,8 kg
4,2 – 6,2 lbs	5,5 – 7,5 lbs	4,2 – 6,2 lbs
Ø75 mm		
FX ₂ – D75 – Up/Down - Low	FX ₂ – D75 – Up - PLUS	FX ₂ – D75 – Up/Down - Through
3,3 – 3,5 kg	4,1 – 4,3 kg	3,3 – 3,5 kg
7,3 – 7,7 lbs	9,0 – 9,5 lbs	7,3 – 7,7 lbs
Ø100 mm		
FX ₂ – D100 – Up/Down - Low	FX ₂ – D100 – Up - PLUS	FX ₂ – D100 – Up/Down - Through
4,8 – 6,9 kg	5,2 – 7,9 kg	4,8 – 5,9 kg
10,6 – 15,2 lbs	11,5 – 17,4 lbs	10,6 – 13,0 lbs

Temperature:
t = max. 70 °C (158 °F)
(extracted air)

Material recovery:
99 % by weight


Vacuum:
p = max. 3500 Pa (0,51 psi)





FX ₂ D50 arms:	FX ₂ D75 arms:	FX ₂ D100 arms:
Recommended		
Q = 50 - 110 m ³ /h	Q = 110 - 240 m ³ /h	Q = 200 - 450 m ³ /h
Q = 29,4 – 64,7 cfm	Q = 64,7 - 140,3 cfm	Q = 117,7 - 264,9 cfm
Maximum		
Q = 142 m ³ /h	Q = 318 m ³ /h	Q = 565 m ³ /h
Q = 83,6 cfm	Q = 187,2 cfm	Q = 332,6 cfm

3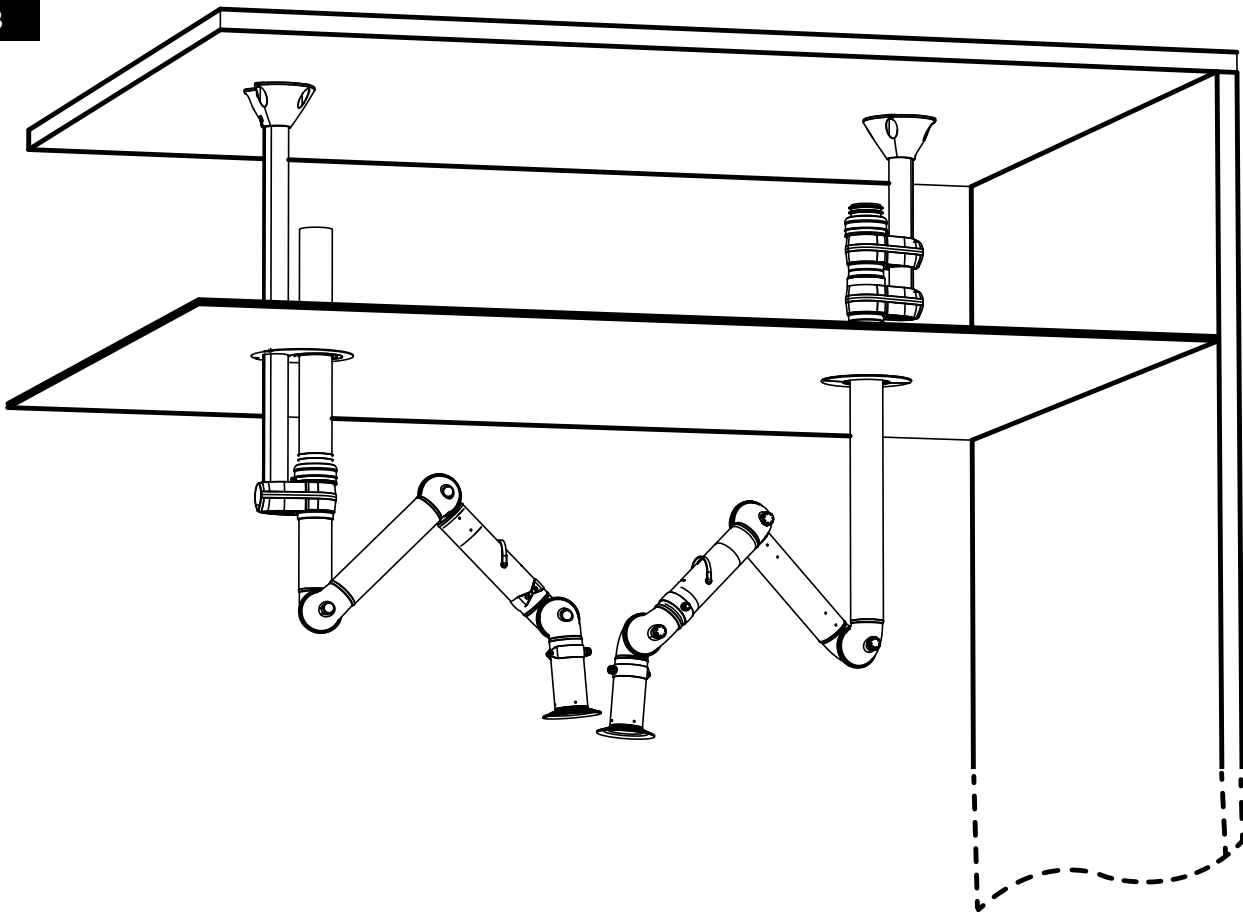


4

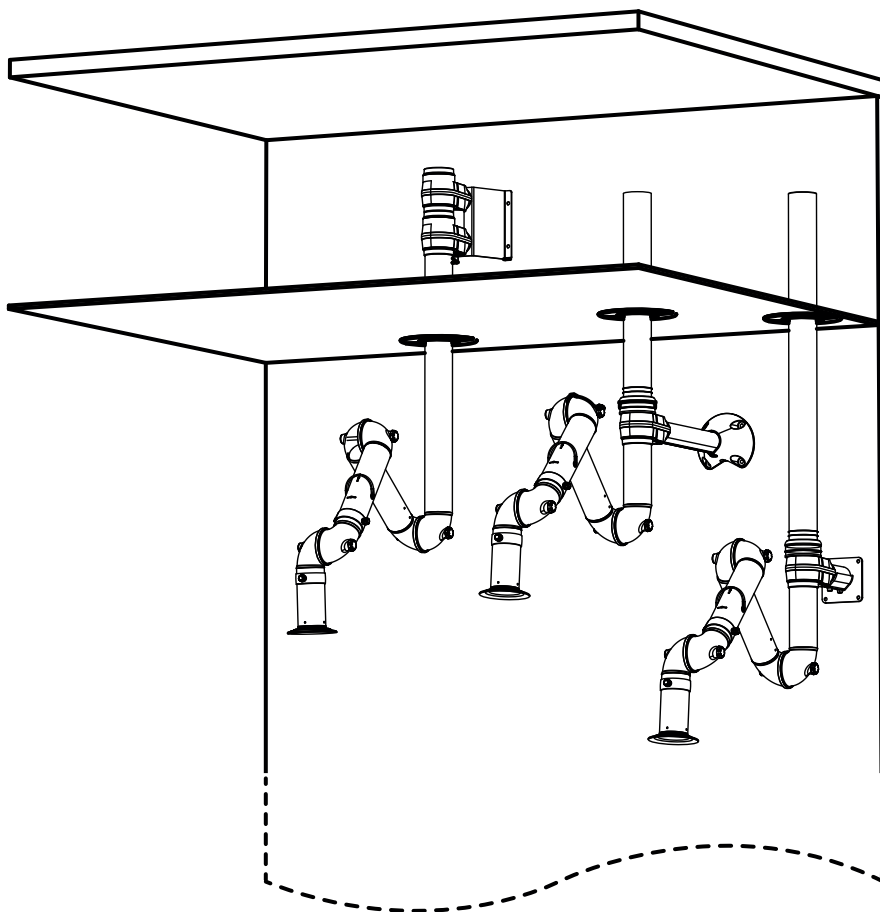
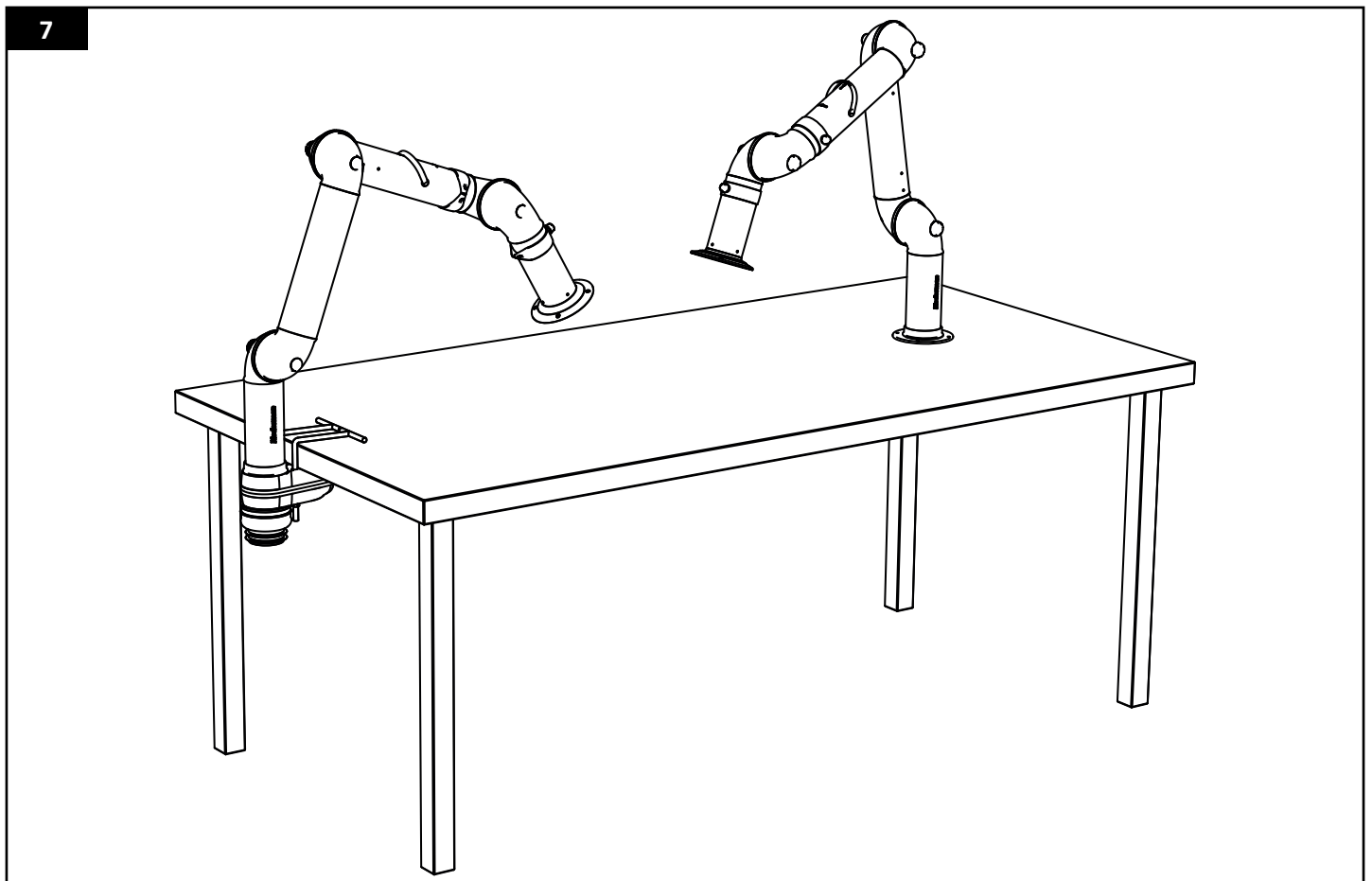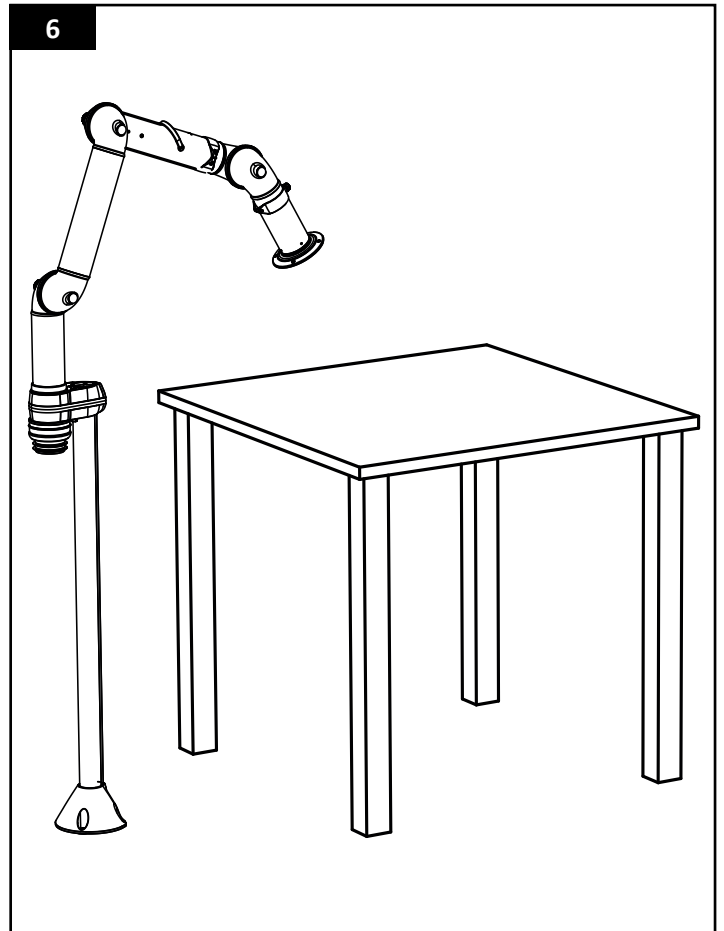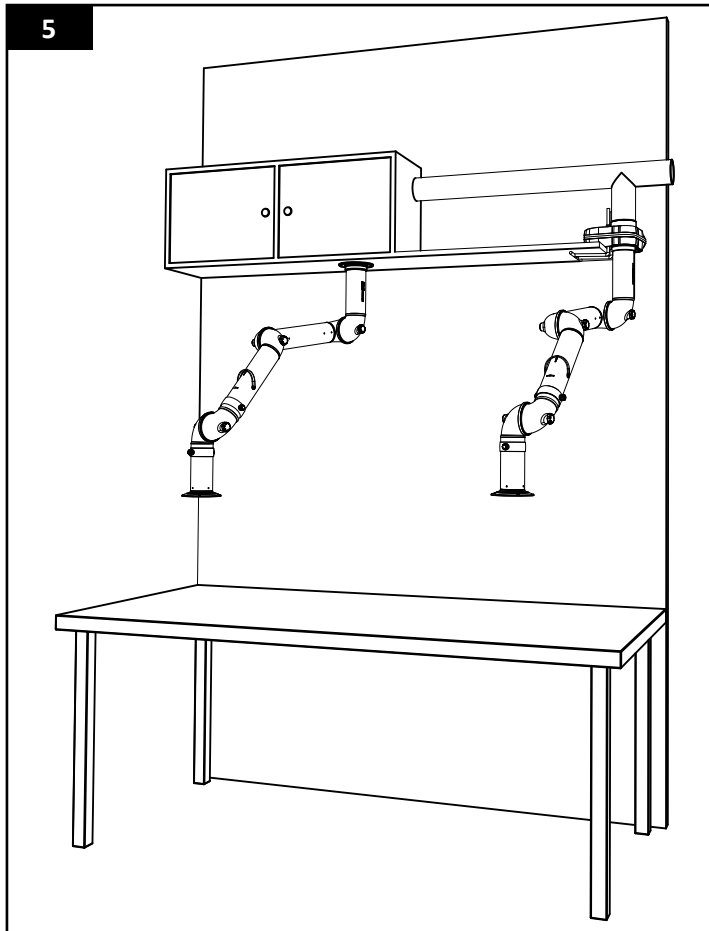
8

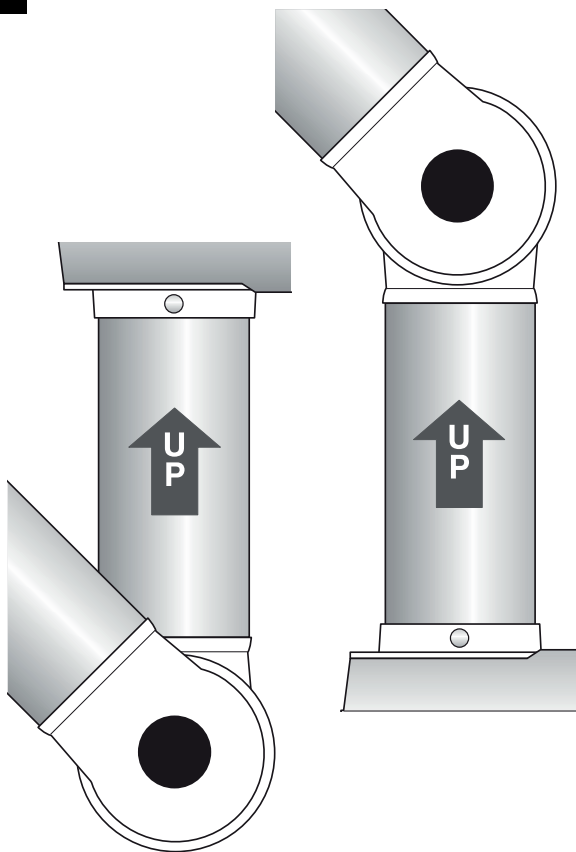9

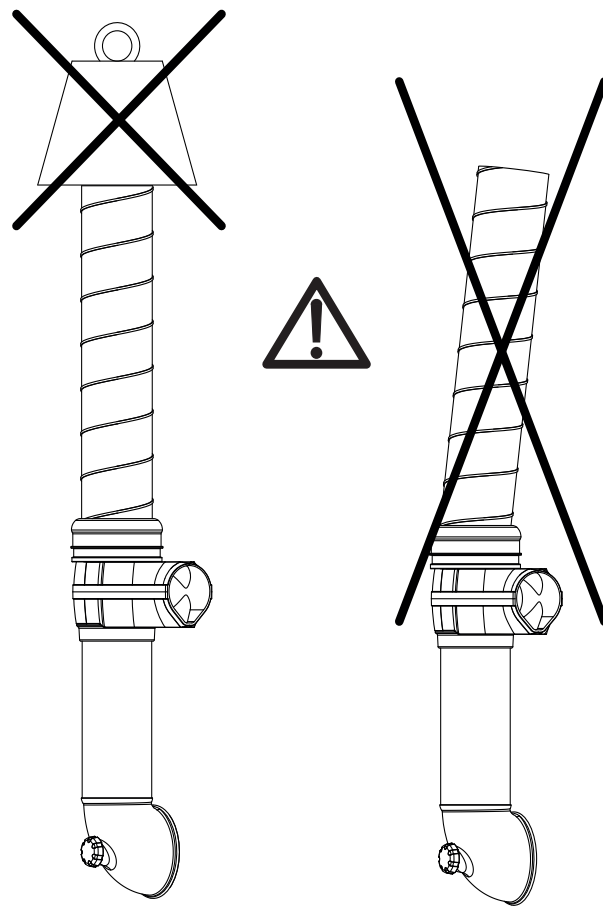
10

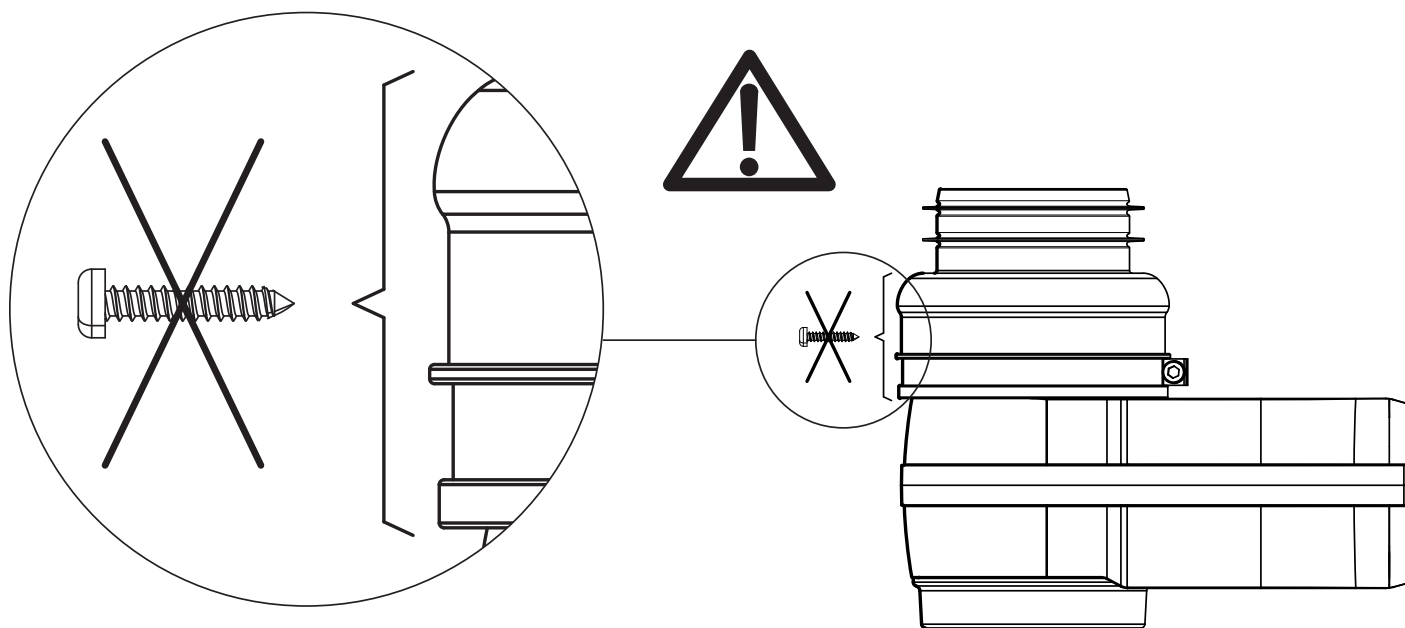
11

SWIVEL AND LINK FUNCTIONS


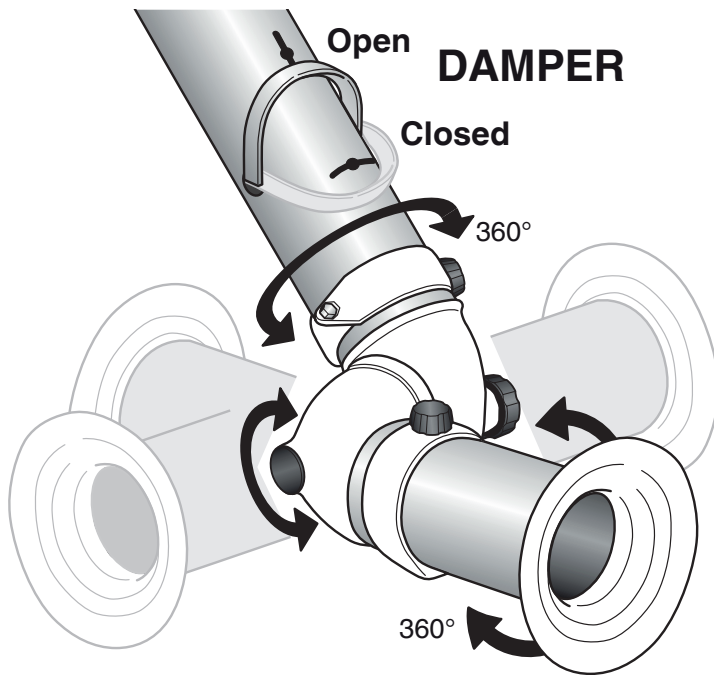



FITTING THE GAS SPRING (FX₂ D100 - L2400)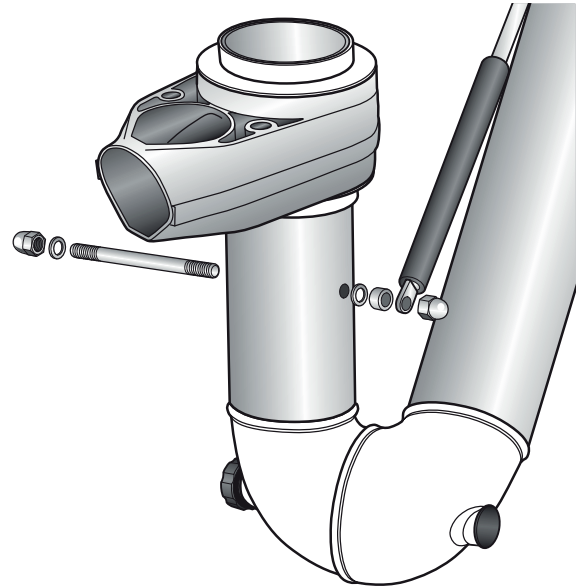

12

i A

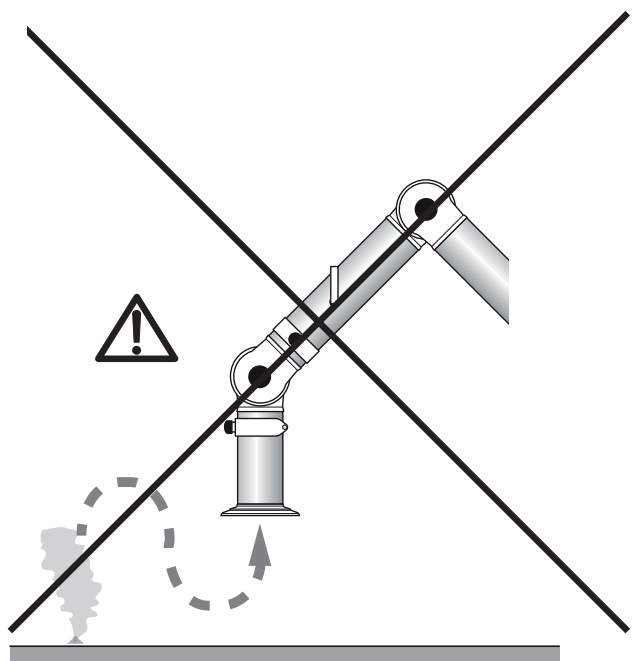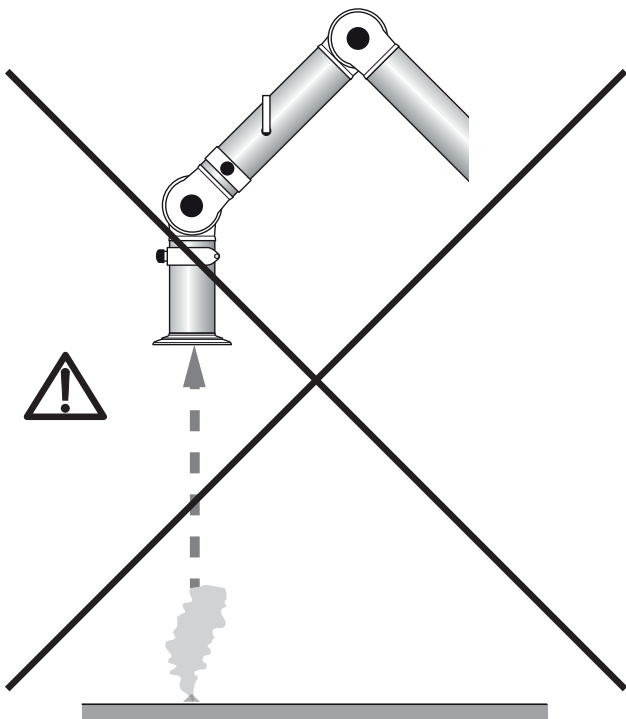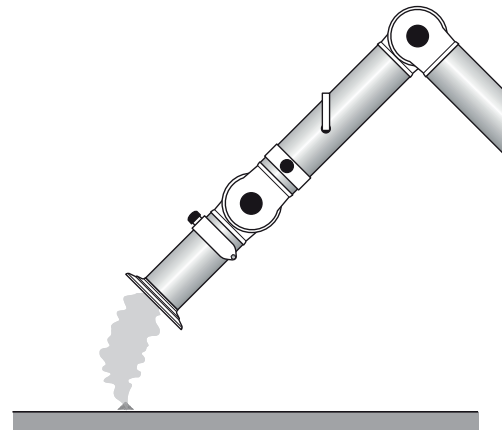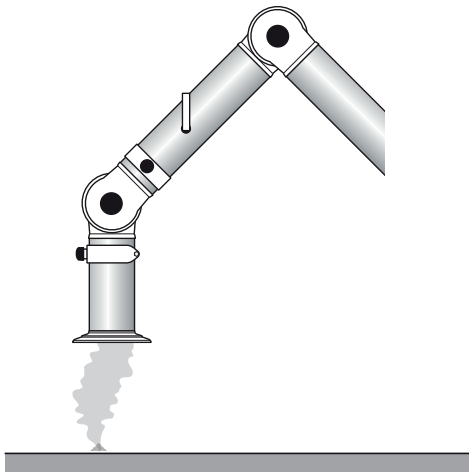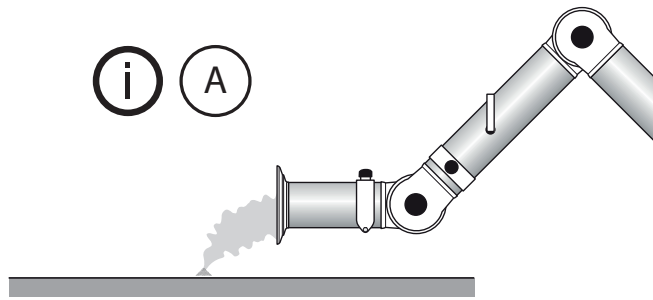English

User Manual

Table of contents

1	Preface.....	14
2	Safety.....	14
3	Description.....	14
3.1	Intended use.....	14
3.2	Technical specifications.....	14
3.2.1	Dimensions.....	14
3.2.2	Technical data.....	14
4	Installation.....	15
4.1	Delivery check.....	15
4.2	Mounting instruction.....	15
4.3	Adjusting the links.....	15
4.4	Damper, swivel and link functions.....	15
5	Operation.....	15
5.1	Arm position.....	15
6	Maintenance.....	15
6.1	Routine inspection and service.....	15
6.2	Spare parts.....	16
6.3	Recycling.....	16
	Appendix A: Installation protocol.....	17
	Appendix B: Maintenance checklist.....	18

1 Preface

Read all product documentation and product identification plate carefully before installation, use, and service of this product. Replace documentation immediately if lost. Nederman reserves the right, without previous notice, to modify and improve its products including documentation.

This product is designed to meet the requirements of relevant EC directives. To maintain this status, all installation, maintenance and repair is to be done by qualified personnel using only Nederman original spare parts and accessories. Contact the nearest authorized distributor or Nederman for advice on technical service and obtaining spare parts. If there are any damaged or missing parts when the product is delivered, notify the carrier and the local Nederman representative immediately.

2 Safety

This document contains important information that is presented as either a warning, caution or note. See the following examples:

WARNING! Risk of personal injury.

Warnings indicate a potential hazard to the health and safety of personnel, and how that hazard may be avoided.

CAUTION! Risk of equipment damage.

Cautions indicate a potential hazard to the product but not to personnel, and how that hazard may be avoided.

NOTE! Notes contain other information that is important for personnel.

3 Description

3.1 Intended use

FX₂ Original: Fumes, vapour

FX₂ Original: Dust

FX₂ Chem: Chemical applications

3.2 Technical specifications

3.2.1 Dimensions

See figure 1.

3.2.2 Technical data

See figure 2.

4 Installation

4.1 Delivery check

If there are any damaged or missing parts when the product is delivered, notify the carrier and the local Nederman representative immediately.

4.2 Mounting instruction

See figures 3, 4, 5, 6 and 7. The extractor arm can be mounted on wall, bench, floor or ceiling. A mounting instruction comes together with the accessories wall/ceiling bracket, hood and bench bracket.

See figure 8. The arm must always be mounted so that the adjusting knobs on the links are on the right hand side.

If an UP-arrow is positioned near the duct connector, the arm must only be installed so that this arrow points upwards.

See figure 9. Make sure the arm and the duct are properly connected.

See figure 10. Do not use a screw with that part of the arm.

4.3 Adjusting the links

The links are factory set for normal working position. Always check and adjust the links so they will suit the users preference.

4.4 Damper, swivel and link functions

See figure 11.

5 Operation

5.1 Arm position

See figure 12. Always place the arm as close as possible to the source. When possible, place the arm sideways to the source for best capture, to avoid interfering with the work process, and to minimize the effect from cross draft. See figure 12(A).

6 Maintenance

CAUTION! Risk of equipment damage.

Use only Nederman original spare parts and accessories.

6.1 Routine inspection and service

Follow the list in. *Appendix B: Maintenance checklist* to routinely inspect, and repair or replace worn and damaged parts on the inside and outside of the product.

6.2 Spare parts

**CAUTION! Risk of equipment damage.**

Use only Nederman original spare parts and accessories.

Contact your nearest authorized distributor or Nederman for advice on technical service or if you require help with spare parts. See also www.nederman.com.

Ordering spare parts

When ordering spare parts always state the following:

- Part number and control number (see the product identification plate).
- Detail number and name of the spare part (see www.nederman.com).
- Quantity of the parts required.

6.3 Recycling

The product has been designed for component materials to be recycled. Its different material types must be handled according to relevant local regulations. Contact the distributor or Nederman if uncertainties arise when scrapping the product at the end of its service life.

Appendix A: Installation protocol

Copy the installation protocol, fill it in and save it as a service record.

For values, note the value in the result column, otherwise a tick will suffice if the item has been performed or considered.

NOTE! If a value is outside its accepted level, or a result is incorrect or missing, fix the problem before initial start-up and normal operation.

Unit No.	Date:	
	Performed by:	

[illegible]

Copy the maintenance checklist, fill it in and save it as a service record.

For values, note the value in the result column, otherwise a tick will suffice if the item has been performed or considered.

NOTE! If an inspection result (for example, a measured value) differs significantly from a previous result, find the cause for the difference.

Unit No.	Date:	
	Operating hours:	
	Performed by:	

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